

Frozen Material Management System

«**Selen Freezer v.1.03**»

Server installation/deinstallation instructions

Mannheim 2009

Installation of Selen Freezer Server

Program installation:

1. Visit the company website under <http://www.selensystems-online.de> or <http://www.selensystems.com>.
2. In the section **Products** choose the configuration of Selen Freezer you are interested in and click on the link **More**.
3. You will be taken to the page with detailed description of the product. On the left hand side below the picture of the product you will find the **Download** link. Click on this link and choose the folder where you want the file to be saved in.
4. After downloading execute **Freezer_Server.exe** will extract installation files (Figure 1).

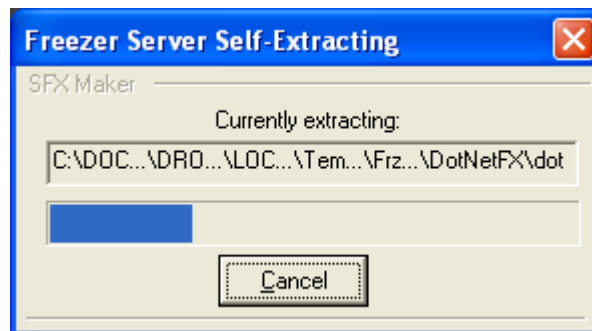


Figure 1. **Selen Freezer Server** file extraction window.
After extraction of all necessary files the installation starts (Fig 2).

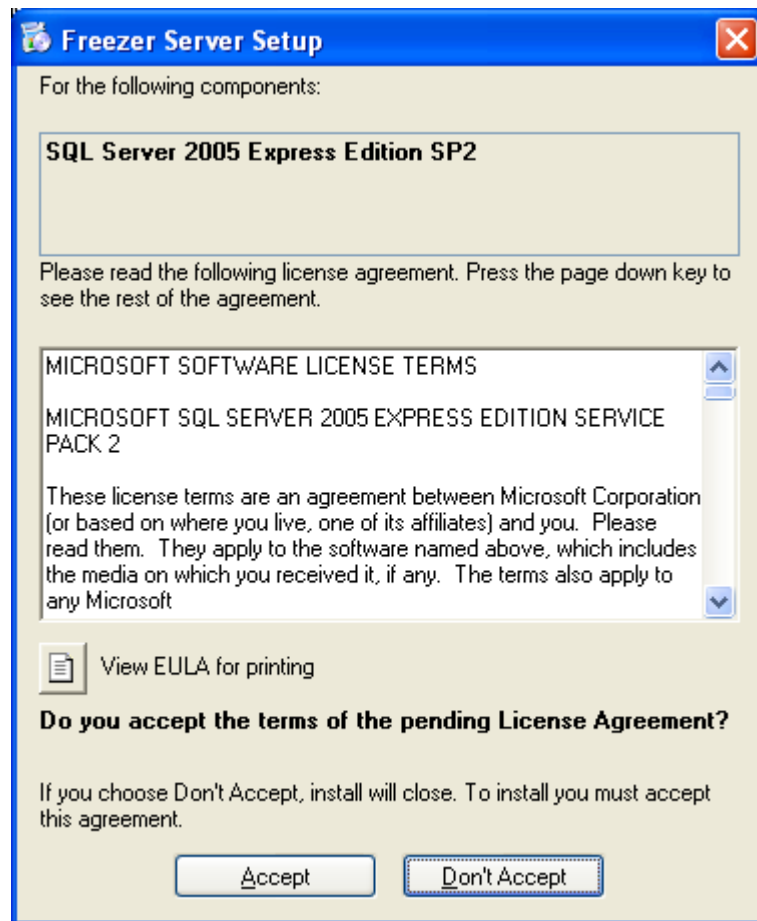



Figure 2. Selen Freezer Server installation window

5. As a first step installation of **SQL Server 2005 Express Edition SP2**.

Important!

Installation of **SQL Server 2005 Express Edition SP2** is necessary for Selen Freezer Server functioning.

Using **View EULA for printing**  you can view the license agreement in **rtf** and print it out.

If you do not want to accept the license agreement click **Don't Accept** button. This will terminate the installation of **SQL Server 2005 Express Edition SP2** and consequently the installation of **Selen Freezer Server**.

Accept the license agreement by pressing **Accept** button. Installation of **SQL Server 2005 Express Edition SP2** will start and the window presented on Figure 3 will appear.

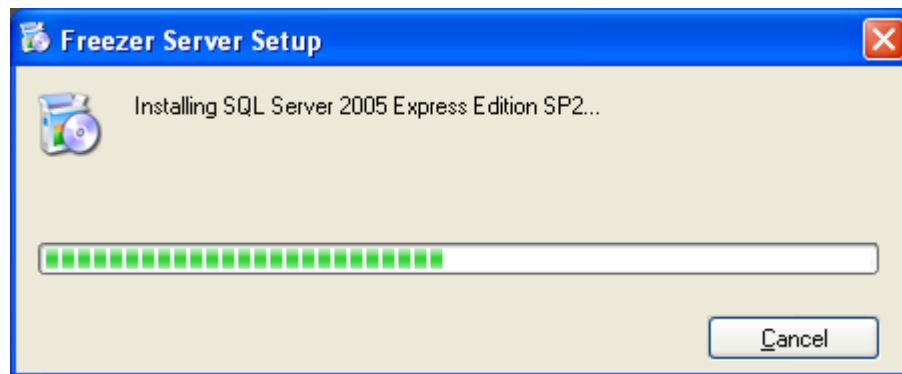


Figure 3. Installation status of **SQL Server 2005 Express Edition SP2**

6. After completion of **SQL Server 2005 Express Edition SP2** c installation you a window presented on the Figure 4 will appear.

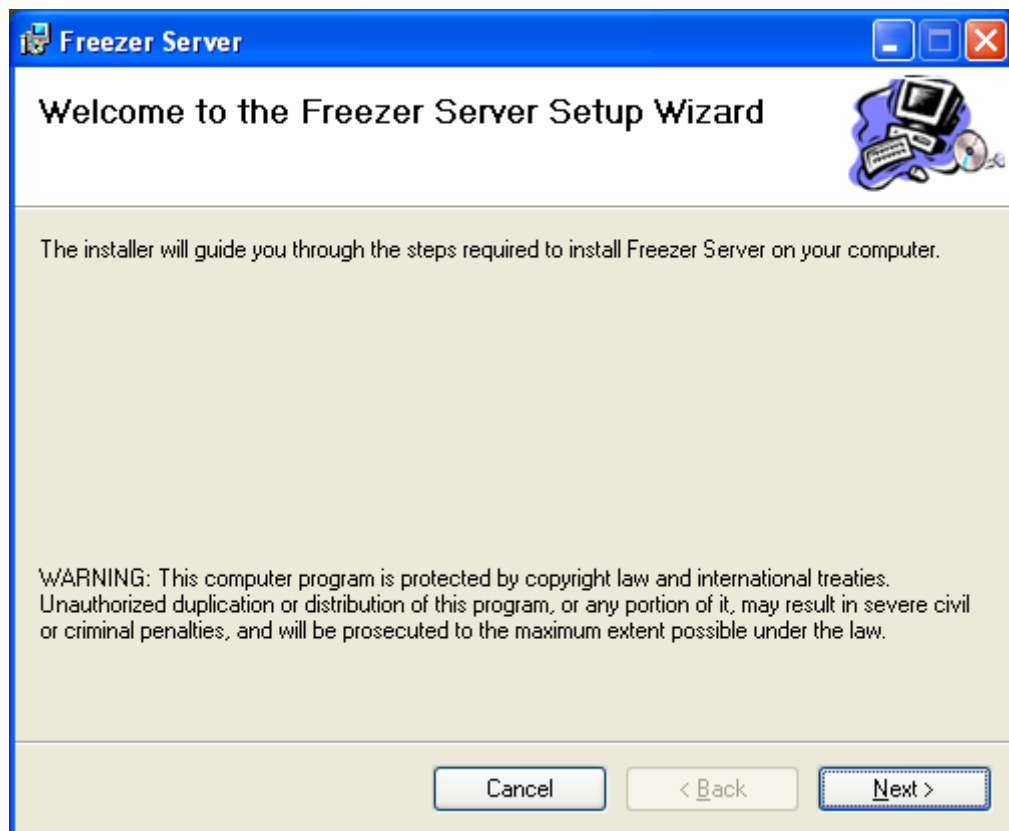


Figure 4. Selen Freezer Server installation window.

7. Press **Next** to proceed to Selen Freezer Server installation. Window presented on figure 5 will appear.

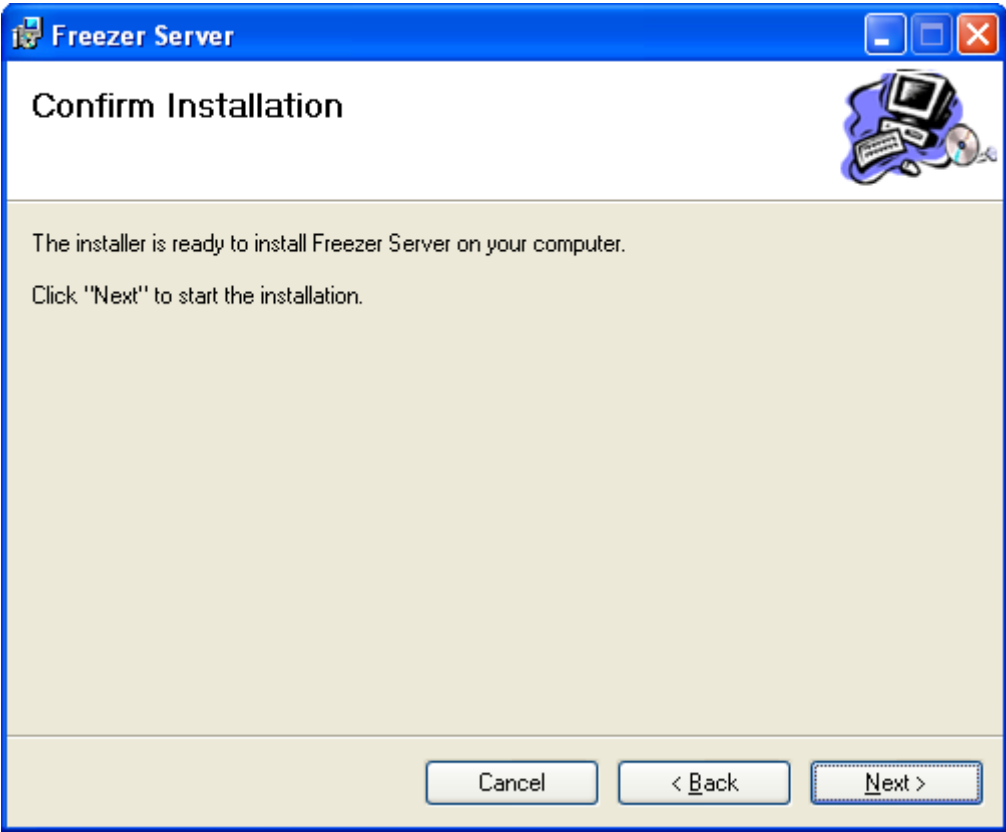


Figure 5. Selen Freezer Server installation window

8. Press **Next** to proceed to Selen Freezer Server installation.

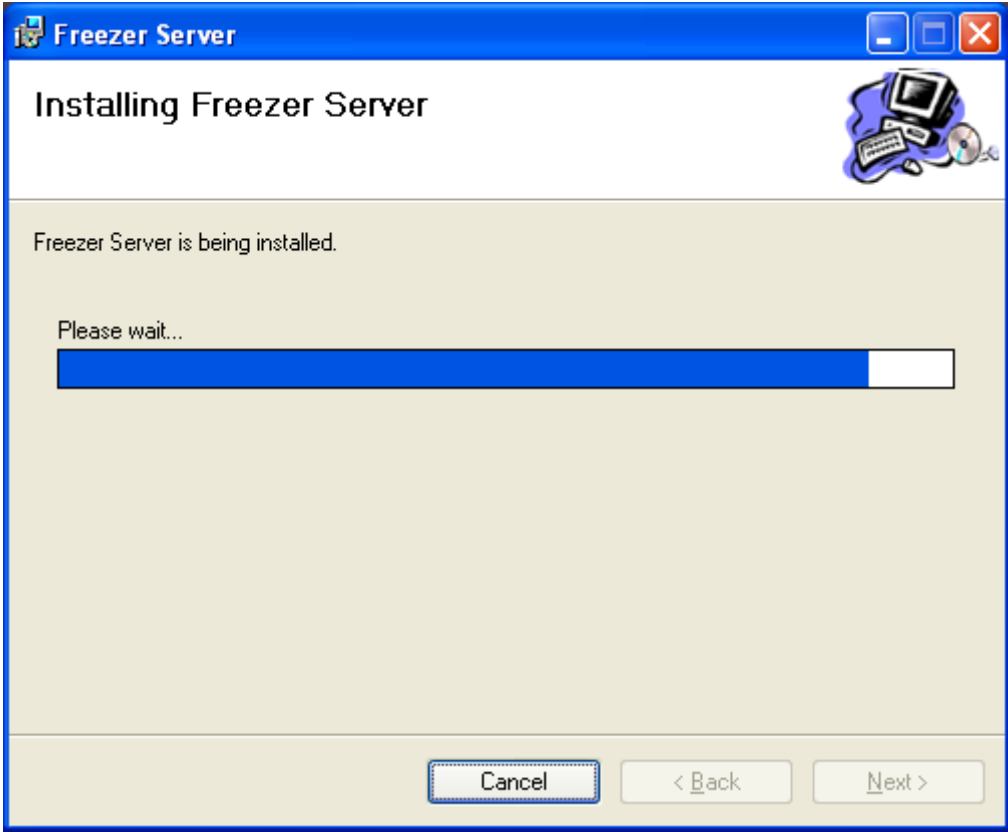


Figure 6. Window showing Selen Freezer Server installation progress.

9. After completing the installation you will see the message, presented on the figure 7.

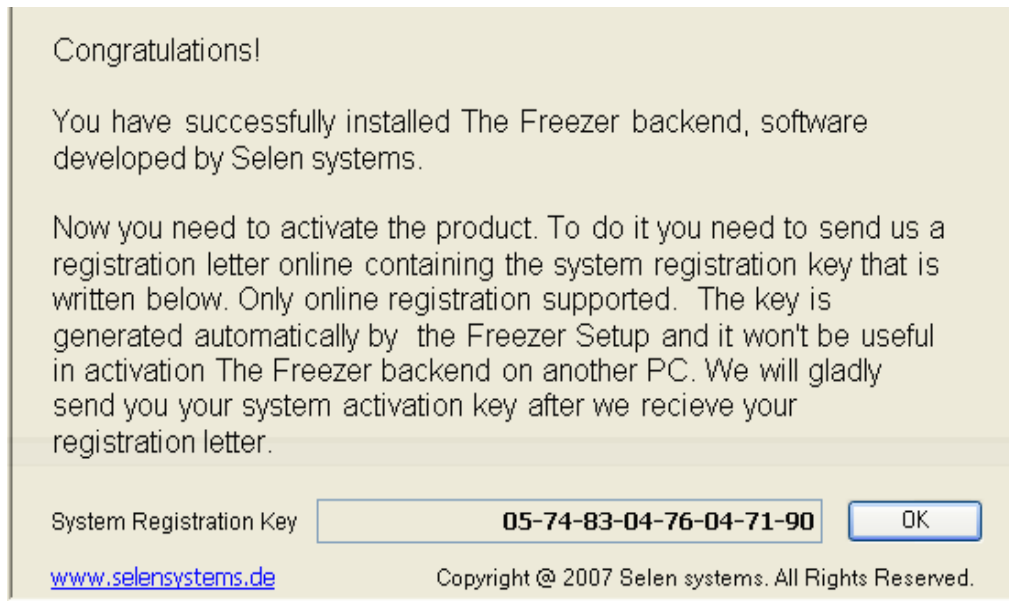


Figure 7. Window with system registration key necessary for Selen Freezer licensing.

Important!

You will need the System Registration Key for obtaining a license file.

Press **OK** button to complete the installation. This will be confirmed by a message presented on the figure 8.

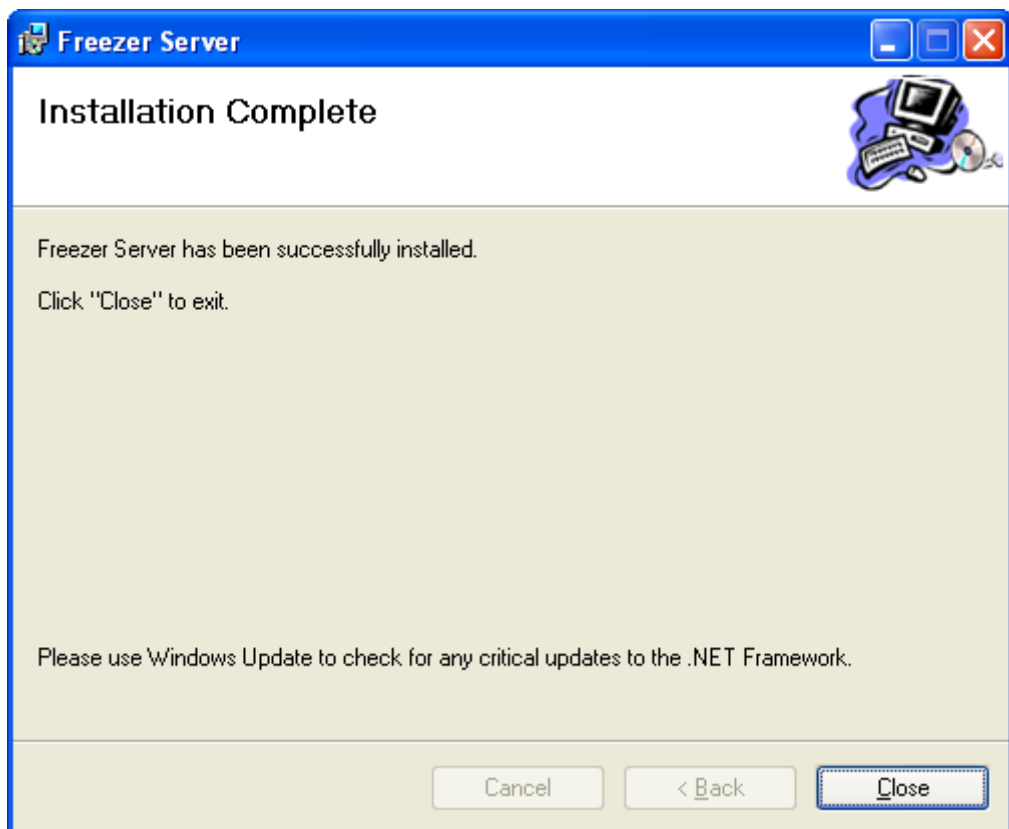


Figure 8. Successful installation confirmation

Press **Close** button to finalize **Selen Freezer Server** installation procedure.

Note

If you want to cancel the installation you can do it before the appearance of System Registration Key by pressing **Cancel** button. If you do so a message presented on figure 9 will appear.

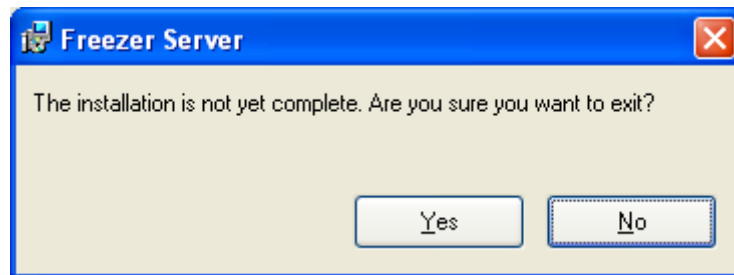


Figure 9. Cancellation of **Selen Freezer Server** installation

If you press **Yes** button the installation will be cancelled and message presented on figure 10 will appear.

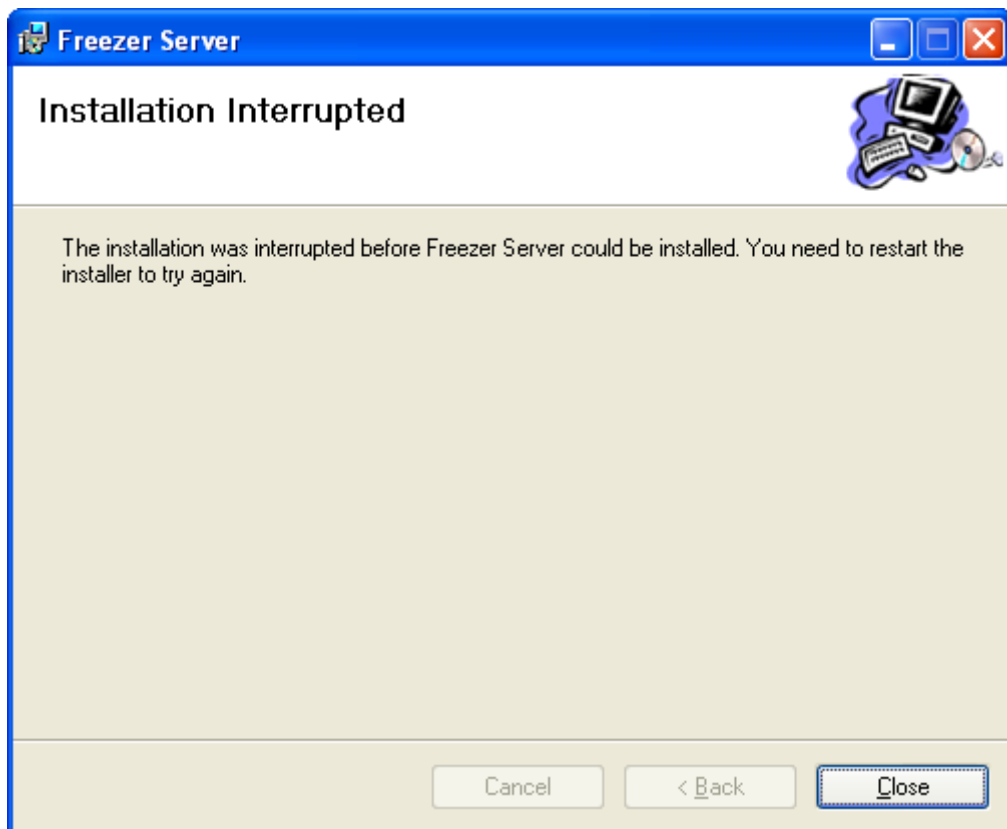


Figure 10. Cancellation of **Selen Freezer Server** installation

Important!

After the installation of **Selen Freezer Server** you can find 3 highly important files in the program folder created during the installation (usually **C:\Program Files\Selen Systems\Freezer Server**). Below you will find the description of these files.

1. File **Key.txt** – is a text file, containing the System Registration key (presented on figure 7) you may need it if you did not save the code during the installation.

2. File **licence.frz** is a demo version license file that has to be copied in the “\Selen Systems\Freezer” folder (folder where the client program is located) if you want to use demo version of the software.

Note

Typical path to client program is **C:\Program Files\Selen Systems\Freezer**. As soon as you copy **licence.frz** into this folder you can use Selen Freezer in demo mode. Demo mode allows generation of 2 users and 2 freezers. To be able to create more freezers or users you have to purchase a license.

3. File **BackUp.exe** is a backup utility that automatically generates backup of **Selen Freezer database**. To test whether backup utility is correctly installed you have to perform following steps:

Important!

Do not remove **BackUp.exe** utility; this will significantly impair the data safety.

Deinstallation of Selen Freezer Server

- Go to start menu;
- Open system settings;
- Choose Software installation/deinstllation;
- Find **Freezer Server**;
- Choose remove option
- Confirm that you want to deinstall the program by pressing Yes button

Note

Please note, that this will remove Selen Freezer Server only! If you do not need **SQL Server 2005 Express Edition SP2**, you will have to remove it separately.